

Multi-Laser Pointing Stability and Side-Lobe Suppression in a Compact Integrated Excitation Module for Flow Cytometry and Microscopy

Luo, N. ¹, He, H. ¹, Xu, M. ¹, Li, J. ¹, Huang, R. ¹, Bergles, E. ¹, Gebhart, D. ¹

¹ Pavilion Integration Corporation



OBJECTIVES & CORE PERFORMANCE

1. Abstract & System Architecture

High-performance fluorescence instrumentation demands perfectly co-aligned, multi-wavelength illumination. Traditional loose-component setups are highly susceptible to thermal misalignment, large structural footprints, and elevated spatial noise.

This work evaluates a compact, integrated multi-laser module unifying **405 nm, 488 nm, 561 nm, and 638 nm** channels into a single, shared optical engine. This design yields wavelength-combined delivery engineered directly for parallel microfluidic flow-cells or advanced microscopy applications while strictly preserving alignment and beam quality.

Co-linear Delivery: Dichroic combiner elements lock alignment mechanically inside a ruggedized housing.

Target Interface: Optimized for tight, reliable delivery into microfluidic streams.

2. Baseline Beam Verification

Initial verification confirmed that the module's focal plane profiles and power attributes easily meet strict engineering specifications over extended operation.



L*W*H: 100*135*50 - 180*155*52.2mm

POINTING STABILITY & PRODUCTION CAPACITY

3. Spatial Alignment Drift Dataset

Spatial alignment drift compromises quantitative precision and causes inter-channel cross-talk in flow cytometry. Beam pointing stability was monitored across 39 discrete operational intervals per channel. Total deviation on both axes was strictly bound to a few micrometers.

	Pointing Stability (um)					
	488nm Channel		638nm Channel		405nm Channel	
	X	Y	X	Y	X	Y
Average \bar{P}	2.24	3.38	1.61	2.94	2.22	3.74
Standard Deviation σ	0.42	0.44	0.46	0.68	0.55	0.83
MAX	3.06	4.31	2.73	4.28	3.45	5.74
Min	1.39	2.31	0.64	1.82	1.30	2.42
$\bar{P} + 3\sigma$	3.49	4.70	2.98	4.98	3.86	6.23
UCL (3 σ + Test System uncertainty)*	4.49	6.20	3.98	6.48	4.86	7.73
Cpk	1.8	3.1	1.57	1.36	1.56	1.46

Statistical Analysis (N=39 samples/channel)

4. Manufacturing Process Capability

To ensure this compact optical design transfers seamlessly to high-volume manufacturing, process capability metrics were derived from the experimental dataset.

PROCESS CAPABILITY INDICES (Cpk PER CHANNEL)

488 nm Channel: [] High (-> 1.0)

638 nm Channel: [] High (-> 1.0)

405 nm Channel: [] High (-> 1.0)

Manufacturing Note: All active optical channels demonstrated exceptionally high process capability indices (Cpk). The qualified probability of units remaining strictly within target boundaries approaches unity, validating production-ready consistency.

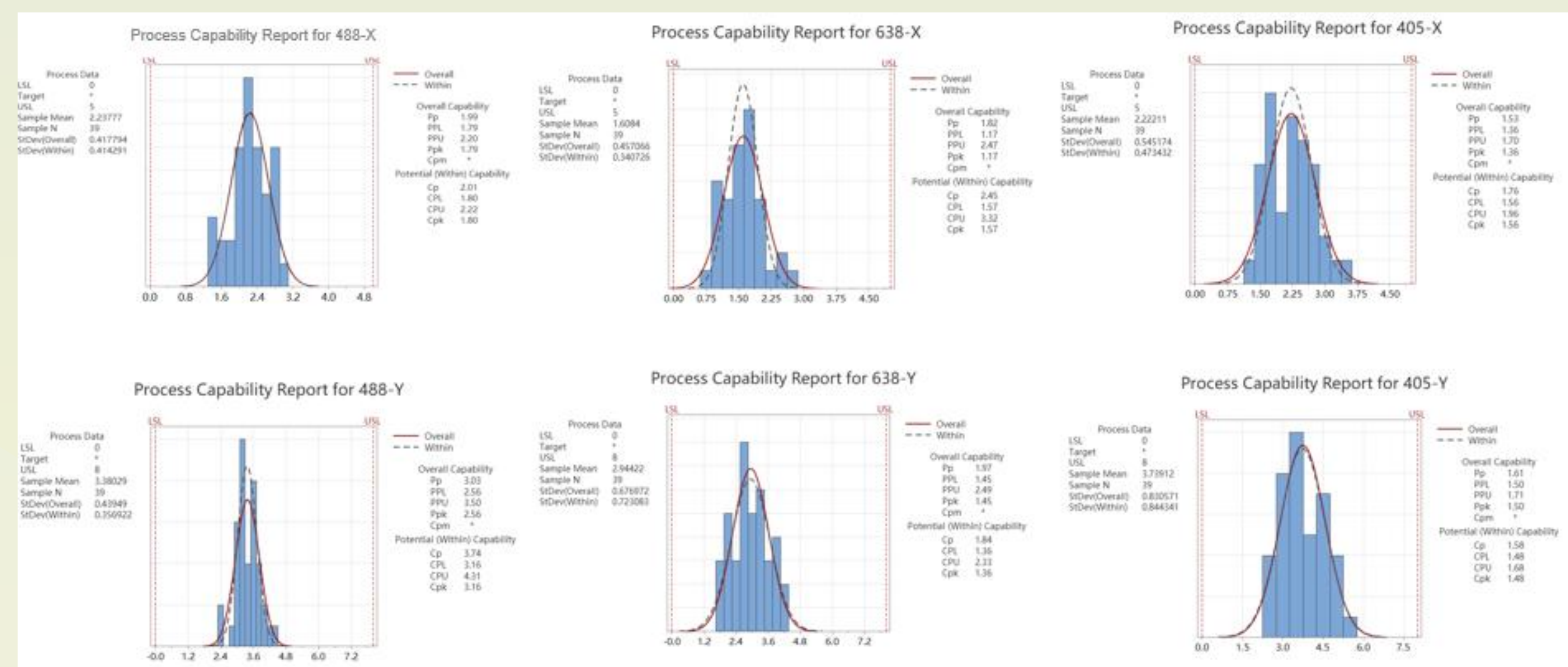


Figure 1: Process Capability of Pointing Stability Data. N=39

SPATIAL NOISE SUPPRESSION & APPLICATION


5. Side-Lobe Suppression Profiles

Spurious spatial side-lobes introduce background noise and stray light artifacts during cellular interrogation. The integrated optical module was benchmarked directly against a traditional loose-component comparator system.

SPATIAL SIDE-LOBE CONTENT COMPARISON (%)

Lower percentage represents a cleaner optical profile

Integrated Module (Typical) | [] < 3%
 Integrated Module (Max Spec) | [] < 5%
 Comparator System (Standard) | [] 10% (Spec Limit)

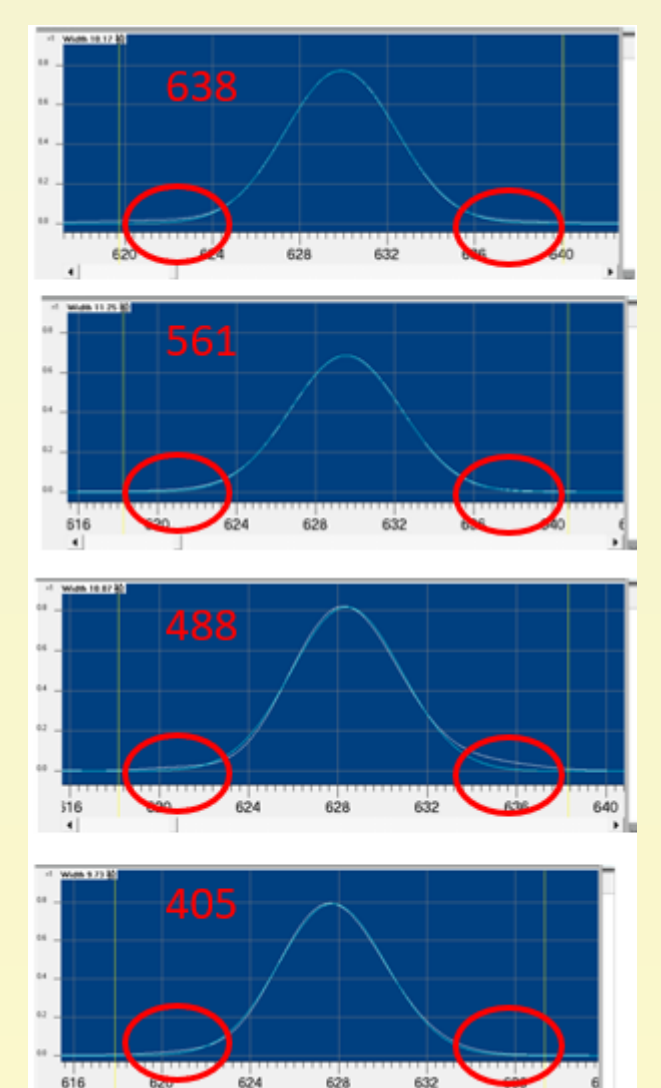

Performance: The integrated module successfully restricts side-lobe content to **below 5%**, and typically **below 3%**.

Advantage: This represents a major engineering improvement over the standard **10%** maximum specification limit typical of unintegrated competitor systems.

CONCLUSIONS

The performance data validates that this integrated multi-laser platform delivers the spatial precision and structural reliability required for demanding fluorescence instrumentation:

- **Sub-Micrometer Control:** Low micro-scale spatial drift guarantees highly repeatable beam targeting.
- **Optically Clean Output:** Side-lobe containment under 3% dramatically lowers the baseline instrumental noise floor.
- **Production-Ready Layout:** High process validation indices (Cpk) verify robust, volume-reproducible assembly.

REFERENCES

1. Clinical & Laboratory Standards Institute (CLSI) H62: Validation of Automated Systems for Flow Cytometry: <https://clsi.org/shop/standards/h62/>
2. ISO9001 & ISO13485 certifications: <https://www.pavilionintegration.com/resources/>
3. ANSI Z136.1 and IEC 60825-1 standards compliance

ACKNOWLEDGMENTS & CONTACT

Pavilion Integration Corporation,
 2528 Qume Drive, STE 1
 San Jose, CA 95131 USA
 Tel. +1(408)453-8801
www.pavilionintegration.com

